

ULTA Prime GF

ULTA™ Prime GF liquid filter cartridges are utilized for the clarification, stabilization and bioburden reduction of aqueous solutions, media and biologicals. These filters have a high dirt holding capacity and exhibit exceptional flow performance compared to polypropylene filters. The hydrophilic nature of ULTA Prime GF filter cartridges also makes them more suitable for gravity fed systems.

ULTA Prime GF utilizes a glass microfiber filter medium encased within an upstream polypropylene mesh and a downstream non-woven filter support material. ULTA Prime GF filter cartridges are dimensionally stable with no media migration. The pleat pack is supported by an inner polypropylene core and outer polypropylene cage, heat bonded to polypropylene end caps.

Features

- Micron rating range from 0.6 to 10 micron
- Wide range of end caps to allow retrofitting of existing systems
- High capacity filter media giving microbial retention
- Thermal bonded construction
- High filtration area



Figure 1. ULTA Prime GF cartridges and capsule.

Materials of construction

Filtration Membrane: Glass Microfiber

Upstream Support: Polypropylene

Downstream Support: Polypropylene

Inner Support Core: Polypropylene

Outer Protection Cage: Polypropylene

End Caps: Polypropylene

Endcap Insert: 316 Stainless Steel

Standard O-rings/gaskets: Silicone/EPDM

Capsule Body: Polypropylene

Capsule Vent Seals: Silicone



Typical applications

- Prefiltration of biopharmaceuticals including tissue culture media and protein solutions
- Ideal for biologicals such as serum and veterinary vaccines
- Provides optimal protection for final membrane filters
- Ideal for batch or small volume processes

Food and biological safety

Materials conform to the relevant requirements of 21CFR Part 177 and current USP Plastics Class VI - 121°C and ISO10993 equivalents. Low concentrations of surfactant may be present.

Effective filtration area

Up to 0.6 m² (6.3 ft²) per 250 mm (10" module).

Retention characteristics

The retention characteristics of ULTA Prime GF have been determined through controlled laboratory tests challenging with a standard aqueous suspension of ACFTD (AC Fine Test Dust) using on-line laser particle counters.

Media Code β ratio	Micron Rating at Various Efficiencies				
	>99.99% 10000	99.98% 50000	99.90% 1000	99% 100	90% 10
.60	0.60	0.57	0.54	0.32	0.20
1.0	1.00	0.95	0.90	0.70	0.50
003	3.00	2.80	1.80	1.00	0.70
005	5.00	4.70	4.50	3.50	1.00
007	7.00	6.70	6.30	4.50	2.50
010	10.00	8.00	7.00	4.80	2.80
020	20.00	16.00	14.00	10.00	6.00

Recommended operating conditions

Up to 70 °C (158 °F) continuous operating temperature and higher short-term temperatures during CIP to the following limits:

Temperature		Maximum forward dP	
°C	°F	(bar)	(psi)
20	68	5.0	72
40	104	4.0	58
60	140	3.0	44
80	176	2.0	29
90	194	1.0	15
>100 (steam)	>212 (steam)	0.3	4

Capsules may be operated up to a temperature of 40 °C (104 °F) at line pressures up to 5.0 bar (72 psig) for liquids and 4.0 bar (58 psi) in air/gas.



Cleaning and sterilization

ULTA Prime GF cartridges can be repeatedly steam sterilized in situ or autoclaved at up to 121 °C (250 °F). They can be sanitized with hot water at up to 90 °C (194 °F) and are compatible with a wide range of chemicals. Capsules can be repeatedly autoclaved up to 121 °C (250 °F).

Recommended rinse volume

Prior to use - 20 liters per 10" cartridge.

Cartridge flow rates

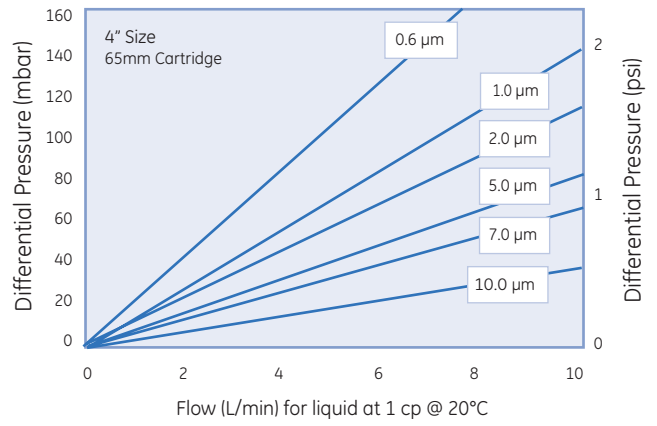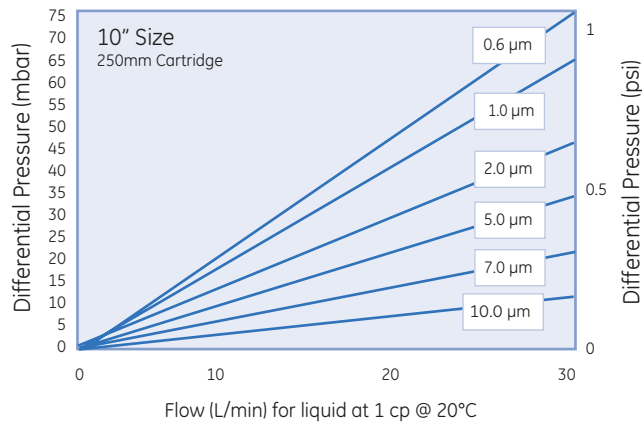

Cartridges

Device Type	Media Type		Special	Detail		Micron Rating		Size	Code Cartridge	O-ring	
Digit 1	2	3	4	5	6	7	8	9	10	11	12
C = Cartridge D = Disc	G	F	Hyphen -	A	Hyphen	9 0 0 0 0 1	6 1 2 5 7 0	H = 5" 1 = 10" 2 = 20" 3 = 30"	0 = Code 0, 222 Double O-ring 7 = Code 7, 226 Locking Tab, Double O-ring/ Fin* (Flat Top 5" Only)	0	E = EPDM S = Silicone

Example: ULTA™ Prime GF, 0.6 µm, 10" 226, Silicone O-ring
CGF-A-96170S

Hyphen	4	7	0	Disc
--------	---	---	---	------

Capsules

Device Type	Media Type		Special	Detail		Micron Rating		Size	Size	End Fittings	
Digit 1	2	3	4	5	6	7	8	9	10	11	12
K = Capsule	G	F	Hyphen -	A	Hyphen	9 0 0 0 0 1	6 1 2 5 7 0	02 = 2" 04 = 4" 06 = 6"		T = 1.5" TC G = SHB* H = 1/2" HB	T = 1.5" TC G = SHB* H = 1/2" HB

Example: ULTA™ Cap PP, 1.0 µm, 4", 1/2 Hosebarb inlet, 1.5 inch TC outlet
KGF-A-0104HT

* Stepped Hose Barb 1/4"

www.gelifesciences.com/filtration

GE Healthcare Bio-Sciences AB
Björkgatan 30, 751 84 Uppsala, Sweden

GE, imagination at work and GE monogram are trademarks of General Electric Company. ULTA is a trademark of GE Healthcare companies.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.

© 2007 General Electric Company - All rights reserved.

GE Healthcare Bio-Sciences AB, a General Electric Company.

GE Healthcare Europe GmbH
Munzinger Strasse 5, D-79111 Freiburg, Germany

GE Healthcare UK Ltd
Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK

GE Healthcare Bio-Sciences Corp
800 Centennial Avenue, P.O. Box 1327
Piscataway, NJ 08855-1327, USA

GE Healthcare Bio-Sciences KK
Sanken Bldg. 3-25-1, Hyakunincho, Shinjuku-ku,
Tokyo 169-0073, Japan

Asia Pacific Tel +65 6275 1830 Fax +65 6275 1829 Australasia Tel + 61 2 9899 0999 Fax +61 2 9899 7511 Austria Tel 01/57606-1619 Fax 01/57606-1627 Belgium Tel 0800 73 888 Fax 02 416 82 06 Canada Tel 800 463 5800 Fax 800 567 1008
Central, East, & South East Europe Tel +43 1 972720 Fax +43 1 97272 2750 Denmark Tel 45 16 2400 Fax 45 16 2424 Finland & Baltics Tel +358 (0)9 512 39 40 Fax +358 (0)9 512 39 439 France Tel 01 69 35 67 00 Fax 01 69 41 96 77
Germany Tel 089 96281 660 Fax 089 96281 620 Greater China Tel +852 2100 6300 Fax +852 2100 6338 Italy Tel 02 27322 1 Fax 02 27302 212 Japan Tel +81 3 5331 9336 Fax +81 3 5331 9370 Latin America Tel +55 11 3933 7300 Fax +55 11 3933 7304
Middle East & Africa Tel +30 210 9600 687 Fax +30 210 9600 693 Netherlands Tel 0800 82 82 82 1 Fax 0800 82 82 82 4 Norway Tel 815 65 555 Fax 815 65 666 Portugal Tel 21 417 7035 Fax 21 417 3184 Russia & other C.I.S. & N.I.S Tel +7 (495) 956 5177
Fax +7 (495) 956 5176 Spain Tel 93 594 49 50 Fax 93 594 49 55 Sweden Tel 018 612 1900 Fax 018 612 1910 Switzerland Tel 0848 8028 12 Fax 0848 8028 13 UK Tel 0800 616928 Fax 0800 616927 USA Tel 800 526 3593 Fax 877 295 8102



imagination at work