

HiPrep 16/10 CM FF HiPrep 16/10 DEAE FF

Instructions

HiPrep™ 16/10 CM FF and HiPrep 16/10 DEAE FF are prepacked, ready to use columns for ion exchange chromatography. They provide fast, preparative separations of proteins and other biomolecules. See below for column characteristics.

Column data

Matrix	6% highly cross-linked spherical agarose	
Mean particle size	90 µm	
Bed volume	20 ml	
Bed height	100 mm	
i.d.	16 mm	
Column composition	Polypropylene	
Recommended flow rate ¹	2–10 ml/min (60–300 cm/h)	
Maximum flow rate ¹	10 ml/min (300 cm/h)	
Maximum pressure over the packed bed during operation, Δp ³	0.15 MPa, 1.5 bar, 22 psi	
HiPrep column hardware pressure limit ³	0.5 MPa, 5 bar, 73 psi	
Storage	+4 °C to +30 °C in 20 % ethanol	
Type of exchanger	CM weak cation	DEAE weak anion
Charged group	-O-CH ₂ COO ⁻	-N ⁺ (C ₂ H ₅) ₂ H
pH stability		
short term	2–14	1–14
working	6–10	2–9
long term	4–13	2–13
Total ionic capacity	0.09–0.13 mmol H ⁺ /ml medium	0.11–0.16 mmol Cl ⁻ /ml medium
Dynamic binding capacity (mg/ml medium) ²		
HSA (M _r 68 000)	N.D.	110
Ribonuclease A (M _r 13 700)	50	N.D.

- 1 Water at room temperature. Flow rate is determined by $v \cdot \eta \leq 10$ ml/min where v =flow rate and η =viscosity.
- 2 Determination of dynamic binding capacity: Samples were applied at 75 cm/h until 50 % breakthrough. Columns: 0.5 x 5 cm. Buffers: 0.05 M Tris, (+2 M NaCl in the elution buffer), pH 7.5 (DEAE), 0.1 M acetate, (+2 M NaCl in the elution buffer), pH 5.0 (CM).
- 3 Many chromatography systems are equipped with pressure gauges to measure the pressure at a particular point in the system, usually just after the pumps. The pressure measured here is the sum of the pre-column pressure, the pressure drop over the medium bed, and the post-column pressure. It is always higher than the pressure drop over the bed alone. We recommend keeping the pressure drop over the bed below 1.5 bar. Setting the upper limit of your pressure gauge to 1.5 bar will ensure the pump shuts down before the medium is overpressured. If necessary, post-column pressure of up to 3.5 bar can be added to the limit without exceeding the column hardware limit. To determine post-column pressure, proceed as follows:

To avoid breaking the column, the post-column pressure must never exceed 3.5 bar.

1. Connect a piece of tubing in place of the column.
2. Run the pump at the maximum flow you intend to use for chromatography. Use a buffer with the same viscosity as you intend to use for chromatography. Note the backpressure as total pressure.
3. Disconnect the tubing and run at the same flow rate used in step 2. Note this backpressure as pre-column pressure.
4. Calculate the post-column pressure as total pressure minus pre-column pressure.

If the post-column pressure is higher than 3.5 bar, take steps to reduce it (shorten tubing, clear clogged tubing, or change flow restrictors) and perform steps 1–4 again until the post-column pressure is below 3.5 bar. When the post-column pressure is satisfactory, add the post-column pressure to 1.5 bar and set this as the upper pressure limit on the chromatography system.

First time use

Ensure an appropriate pressure limit has been set. Equilibrate the column for first time use or after long-term storage by running:

1. 100 ml start buffer (low ionic strength) at 5 ml/min at room temperature (see section "Choice of buffer" for buffer recommendations).
2. 100 ml elution buffer at 5 ml/min at room temperature.
3. 100 ml start buffer at 5 ml/min at room temperature.

These HiPrep columns can be used directly in ÄKTAdesign™ systems without the need for any extra connectors.


Try these conditions first

Flow rate: 5 ml/min at room temperature
Start buffer: See section "Choice of buffer"
Elution buffer: Start buffer + 1 M NaCl
Gradient: 0–100% elution buffer in 200 ml (10 CV)

Equilibration before a new run

Proceed according to steps 2 and 3 in the section "First time use". Extended equilibration may be needed if detergents were included in the eluent.

Please read the back of these instructions for information on optimizing a separation.

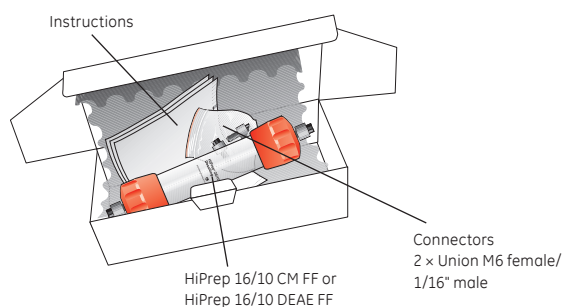

Buffers and solvent resistance

De-gas and filter all solutions through a 0.45 µm filter to increase column life-time.



Daily use

All commonly used aqueous buffers (see "Column data" for recommended pH)
Guanidine hydrochloride, up to 6 M
Urea, up to 8 M



Cleaning

Sodium hydroxide, up to 1 M
Ethanol, up to 70 %
Acetic acid, up to 1 M
Isopropanol, up to 30 %



Avoid

Oxidizing agents
Cationic detergents and buffers (CM)
Anionic detergents and buffers (DEAE)
Phenol

Sample recommendations

Net charge of protein: Positive (CM), negative (DEAE)
Recommended sample load: Not more than 10–20 % of the dynamic capacity (see section "Column data").
Preparation: Dissolve the sample in start buffer, filter through 0.45 µm or centrifuge at 10 000 x g for 10 min



Delivery/storage

The column is supplied in 20% ethanol. If the column is to be stored for more than two days after use, clean the column according to the procedure described under "Cleaning-in-place (CIP)". Then equilibrate with at least 100 ml of 20 % ethanol at a flow rate of 5 ml/min at room temperature.

DO NOT OPEN THE COLUMN!

Choice of buffer

To avoid local disturbances in pH caused by buffering ions participating in the ion exchange process, select a buffer with buffering ions of the same charge as the substituent groups on the ion exchanger.

The start buffer pH should be chosen so that substances to be bound to the ion exchanger are charged, that is, at least 1 pH unit above the isoelectric point for anion exchangers or at least 1 pH unit below the isoelectric point for cation exchangers. Figure 1 and Figure 2 list a selection of standard aqueous buffers.

Table 1 lists suggested volatile buffers used in cases where the purified substance has to be freeze-dried.

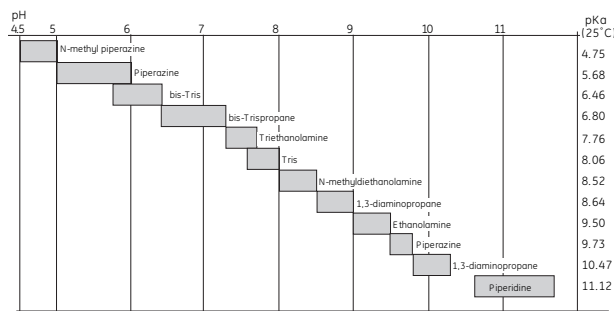


Figure 1. Recommended buffer substances for anion exchange chromatography.

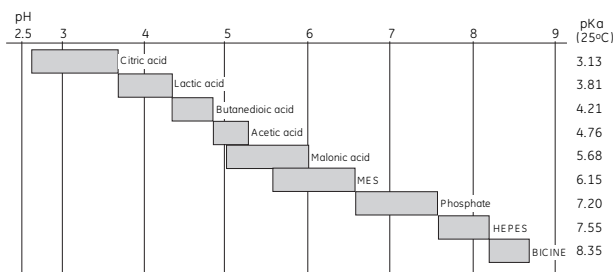


Figure 2. Recommended buffer substances for cation exchange chromatography.

Table 1. Volatile buffer systems.

pH	Substances
2.3–3.5	Pyridine/formic acid
3.0–5.0	Trimethylamine/formic acid
4.0–6.0	Trimethylamine/acetic acid
6.8–8.8	Trimethylamine/HCl
7.0–8.5	Ammonia/formic acid
8.5–10.0	Ammonia/acetic acid
7.0–12.0	Trimethylamine/CO ₂
8.0–9.5	Ammonium carbonate/ammonia
8.5–10.5	Ethanolamine/HCl

Optimization

Perform your first run according to "Try these conditions first". If the obtained results are unsatisfactory, consider the following:

Action	Effect
Change pH/buffer salt (see Figs.1 and 2 for buffers)	Selectivity change,weaker/stronger binding
Change salt, counter ions, and/or co-ions	Selectivity change
Smaller sample loading	Improved resolution
Lower flow rate	Improved resolution
Shallower gradient	Improved resolution, but broader peaks and decreased concentration in fractions

For more information, please check www.gehealthcare.com/protein-purification or refer to the handbook "Ion Exchange Chromatography and Chromatofocusing, Principles and Methods", see ordering information.

www.gehealthcare.com/protein-purification
www.gehealthcare.com

GE Healthcare Bio-Sciences AB
 Björkgatan 30
 751 84 Uppsala
 Sweden



Cleaning-in-Place (CIP)

Regular cleaning

Wash the column with 40 ml of 2 M NaCl at a flow rate of 5 ml/min at room temperature after each run to elute material still bound to the column.

If detergents have been used, rinse the column with 100 ml distilled water followed by 40 ml of 2 M NaCl at a flow rate of 5 ml/min at room temperature.

Re-equilibrate the column with at least 100 ml start buffer at a flow rate of 5 ml/min at room temperature, until the UV baseline and pH/conductivity values are stable.

More rigorous cleaning

Reverse the flow direction and run the following sequence of solutions at a flow rate of 5 ml/min at room temperature.

- 80 ml of a 2 M NaCl solution (removes ionically bound proteins from the column) followed by 50 ml distilled water.
- 80 ml of a 1 M NaOH solution (removes precipitated proteins, hydrophobically bound proteins, and lipoproteins from the column) followed by 50 ml distilled water.
- 80 ml of 70 % ethanol or 30 % isopropanol (removes proteins, lipoproteins, and lipids that are strongly hydrophobically bound to the column) followed by 60 ml distilled water
 or
 30 ml 0.5 % non-ionic detergent in acidic solution (e.g. 0.1 M acetic acid) followed by 100 ml 70% ethanol (to remove the detergent) and 60 ml distilled water.

After cleaning, equilibrate the column before use with approximately 100 ml start buffer at a flow rate of 5 ml/min at room temperature in the normal flow direction.

DO NOT OPEN THE COLUMN!


Trouble shooting

Symptom	Remedy
Increased backpressure over the column	Reverse the flow direction and pump 100 ml elution buffer at a flow rate of 5 ml/min through the column. Return to normal flow direction and run 100 ml buffer at a flow rate of 5 ml/min through the column. If back-pressure is not decreased reverse the flow direction again and follow the rigorous cleaning instructions.
Loss of resolution and/or decreased sample recovery	Clean the column according to the procedure described in the section "More rigorous cleaning".
Air in the column	Reverse the flow direction and pump 100 ml of well de-gassed start buffer through the column at a flow rate of 5 ml/min.

Column performance control


We recommend checking the column performance at regular intervals. Figure 3 describes how to check the function of HiPrep 16/10 CM FF and HiPrep 16/10 DEAE FF.

a) HiPrep 16/10 CM FF

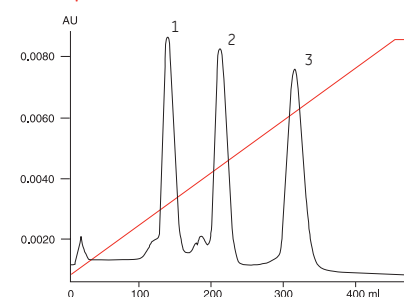

Sample:
 1. Ribonuclease A 1.5 mg/ml
 2. Cytochrome C 0.4 mg/ml
 3. Lysozyme 0.4 mg/ml
 Sample volume: 100 µl
 Gradient: 0–100% elution buffer in 400 ml (20 CV)
 Start buffer: 20 mM Sodium phosphate, pH 6.8
 Elution buffer: 20 mM Sodium phosphate, 0.5 M NaCl, pH 6.8
 Flow rate: 4.0 ml/min, room temperature

b) HiPrep 16/10 DEAE FF

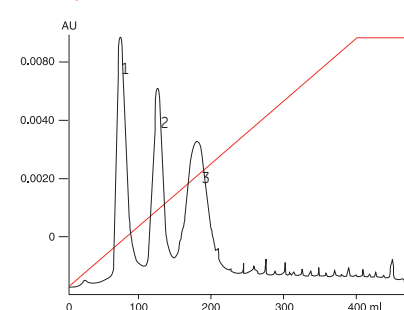

Sample:
 1. α-amylase 2.0 mg/ml
 2. α-Lactalbumin 1.0 mg/ml
 3. Trypsin inhibitor 3.0 mg/ml
 Sample volume: 100 µl
 Gradient: 0–100% elution buffer in 400 ml (20 CV)
 Start buffer: 20 mM Tris, pH 7.5
 Elution buffer: 20 mM Tris, 1.0 M NaCl, pH 7.5
 Flow rate: 4.0 ml/min, room temperature

Figure 3. Typical chromatogram from a function test of (a) HiPrep 16/10 CM FF and (b) HiPrep 16/10 DEAE FF.

Ordering information

Product	No. per pack	Code No.
HiPrep 16/10 CM FF	1 (20 ml)	17-5091-01
HiPrep 16/10 DEAE FF	1 (20 ml)	17-5090-01

Companion products	No. per pack	Code No.
HiPrep 26/10 Desalting	1 (53 ml)	17-5087-01
HiPrep 26/10 Desalting Kit	4 (53 ml)	17-5087-02
HiTrap™ IEX Selection Kit*	7 × 1 ml	17-6002-33
HiTrap DEAE FF	5 × 1 ml	17-5055-01
HiTrap DEAE FF	5 × 5 ml	17-5154-01
HiTrap CM FF	5 × 1 ml	17-5056-01
HiTrap CM FF	5 × 5 ml	17-5155-01

* Contains 7 different ion exchange media

Accessories	No. per pack	Code No.
HiTrap/HiPrep 1/16" male connector to ÄKTAdesign	8	28-4010-81
To connect columns with 1/16" connections to FPLC™ System: Union M6 female/1/16" male*	5	18-3858-01

Literature	No. per pack	Code No.
Handbook, Ion Exchange Chromatography and Chromatofocusing, Principles & Methods	1	11-0004-21
Ion Exchange Chromatography, Media and Column Guide	1	18-1127-31

* Two units (in red polypropylene) are included in the HiPrep package.

GE Healthcare Europe GmbH
 Munzinger Strasse 5, D-79111 Freiburg, Germany

GE Healthcare UK Ltd
 Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK

GE Healthcare Bio-Sciences Corp
 800 Centennial Avenue, P.O. Box 1327, Piscataway, NJ 08855-1327, USA

GE Healthcare Bio-Sciences KK
 Sanken Bldg, 3-25-1, Hyakunincho, Shinjuku-ku, Tokyo 169-0073, Japan

HiPrep, HiTrap, ÄKTAdesign, FPLC, and Drop Design are trademarks of GE Healthcare Ltd. GE tagline and GE Monogram are trademarks of General Electric Company.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. General Electric Company reserves the right, subject to any regulatory and contractual approval, if required, to make changes in specifications and features shown herein, or discontinue the product described at any time without notice or obligation.

© 2006 General Electric Company – All rights reserved.

GE Healthcare Bio-Sciences AB, a General Electric Company.