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The screenshot shows the GE Healthcare Life Sciences login page. At the top left is the GE logo, followed by the text 'GE Healthcare Life Sciences'. Below this is the heading 'Login'. There are two radio button options: 'With User Name' and 'With Activation ID'. The 'With Activation ID' option is selected and highlighted with a red rectangular box. Below the radio buttons is a text input field labeled 'Activation ID'. At the bottom left is a 'Register' link, and at the bottom right is a 'Log in' button.

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This screenshot shows the same login page as the previous one, but now the 'Activation ID' field is filled with the text 'db31-e86f-46d0-413d-b6e8-80b3-c153-999d'. This text is highlighted with a red rectangular box. The 'Log in' button at the bottom right is also highlighted with a red rectangular box. The 'With Activation ID' radio button remains selected.

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The screenshot shows the GE Healthcare Life Sciences License & Delivery Portal. The page has a blue header with the GE logo and navigation links: Home, License Activation, License Support, and Downloads. The main content area is titled 'License & Delivery Portal' and includes a user ID 'db31-e86f-46d0-413d-b6e8-80b3-c153-081d'. There are four main sections: 'Recent Entitlements' with a table of activation IDs and products; 'Recent Files' with a table of file descriptions and dates; 'Announcements' with a test widget; and 'QuickLinks' with links to manage entitlements/licenses and general information/FAQs. The 'Your Downloads' section is highlighted with a red box and contains a download icon and links for 'ALL' and 'UNICORN'.

Activation ID	Product	Last modified
db31-e86f-46d0-413d-b6e8-80b3-c153-081d	UNICORN 7	Jul 27, 2016

Description	Date
UNICORN 7.0.2 Infoletter 29207541AA	Jul 12, 2016
UNICORN 7.0 ZIP	Jun 29, 2016

QuickLinks

- Manage Entitlements
- Manage Licenses
- General Information
- Frequently Asked Questions

© 2015 General Electric Company doing business as GE Healthcare

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	File Description	File Size	File Added	File Name
<input type="checkbox"/>	UNICORN 7.0 ZIP	4.84MB	Jun 29, 2016	↓ UNICORN 7.0.zip

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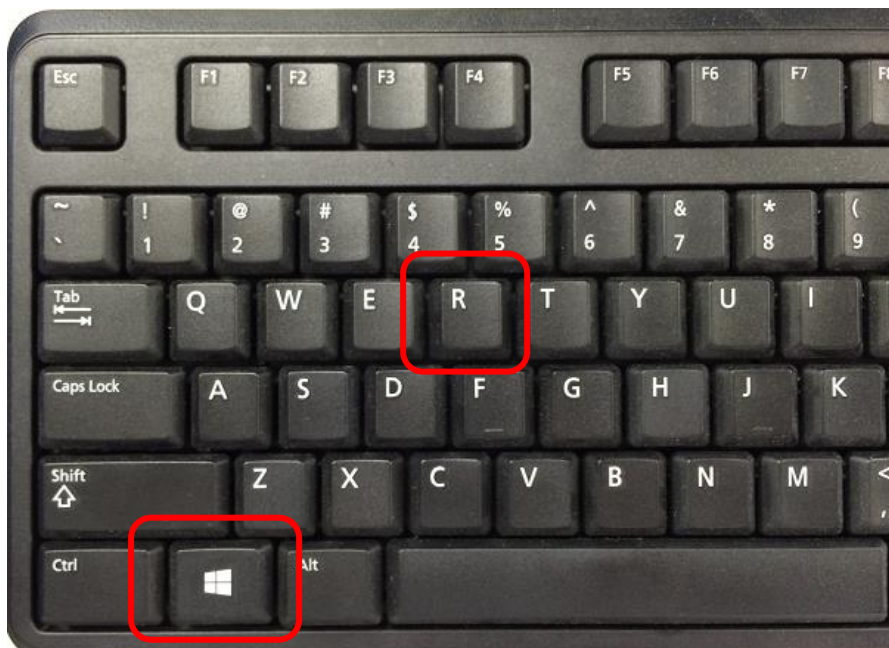
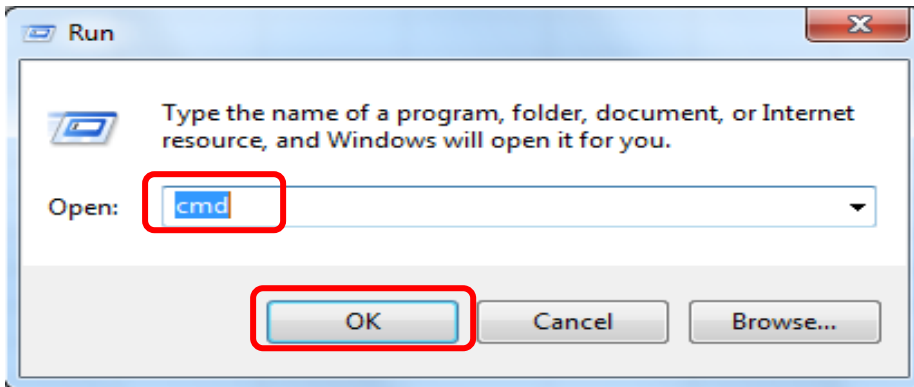
Activate Licenses

db31-e86f-46d0-413d-b6e8-80b3-c153-081d ▼

Please add or select at least one host and click Next to configure counts.

Generate Licenses

Sold to:

Ship to email:

Ship to address:

Selected line items:

Activation ID	Product	Expiration date
db31-e86f-46d0-413d-b6e8-80b3-c153-081d	UNICORN 7	PERMANENT

License Model Attributes

License model type: Nodelocked Uncounted

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Configure Hosts

db31-e86f-46d0-413d-b6e8-80b3-c153-081d ▼

Please enter values for the hosts and click Verify to validate the license. Or click Generate to generate the license.

Configure Hosts


Server hosts: * (no license hosts selected)

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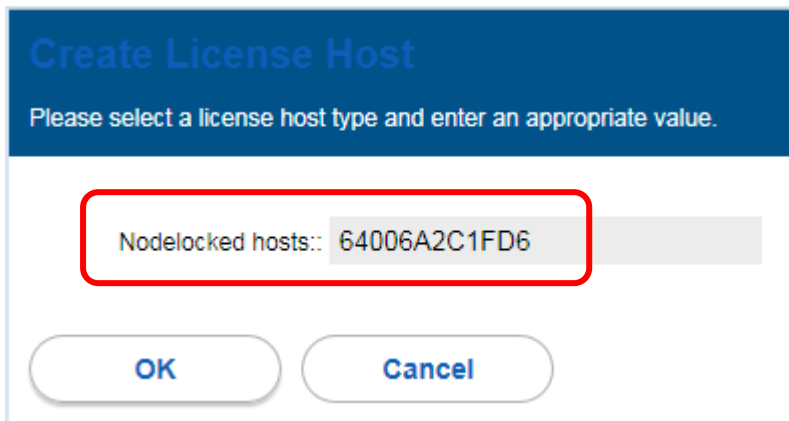
Create License Host
Please select a license host type and enter an appropriate value.

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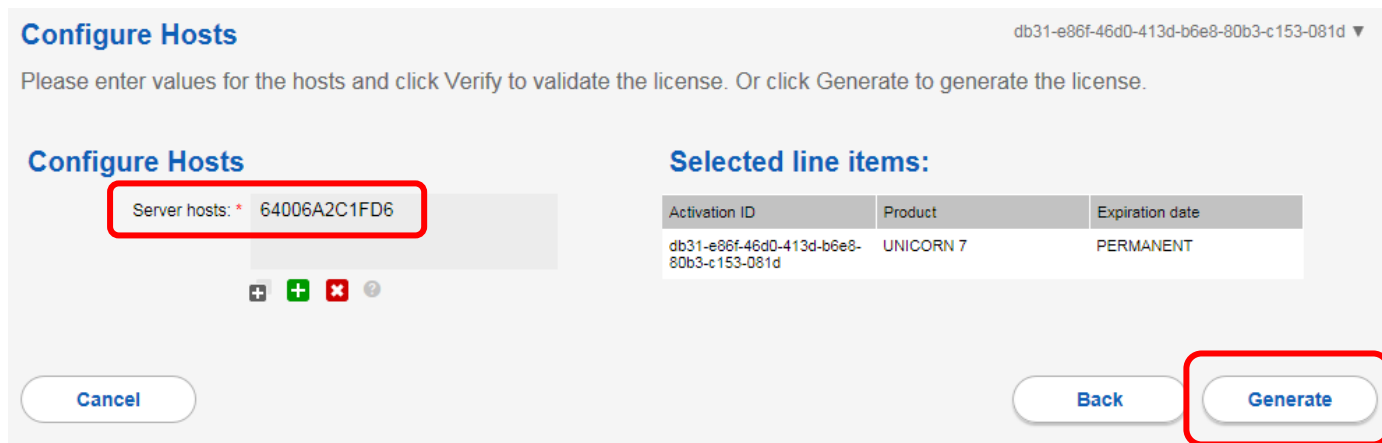Configure Counts 1234-5648-9abc-defg-hijk-lmno-pqrs-tuvw ▼

Please enter values for the counts and click Next to review.

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License Summary

Please review the licenses below and click Complete to return to Manage Entitlements page.

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